

Undercounter Freestanding Freezers (-30°C and -40°C Operation) Temperature Setting



- 1. Press the **SET** button and the **V** button until the **H** parameter is showing. Let go of both buttons.
- 2. Press the button twice. This will display the setting. Press the ET button and the value will be displayed. Change that number using the button to read -10°C then press the ET button. The value will flash on screen and take you to the next setting. At the next setting, continue pressing the button until you reach the HLU setting.
- 3. Once you are at the FLU setting, press the SET button and use the buttons to change the setting to read -15°C. Press the SET button to lock in the setting. The next setting that will appear will be FLL. Press the SET button and using the buttons, change the setting to read -25°C. Press the SET button to lock in the value. Wait 10 seconds for the controller to time out and display the temperature again.
- 4. Once the unit is displaying the temperature, press and hold only the **SET** button until the **L** symbol is flashing. Then use the **SET** buttons to set your unit to **-22°C**. Press the **SET** button again to lock in the temperature value. You are now finished.